

Unpacking the P-737.xSL

NOTICE



Impermissibly high forces on the motion platform!

Impermissibly high forces that are applied to the motion platform can damage the stage.

- Avoid forces on the motion platform of the stage.
 - Only hold the stage externally by the base body.
-

1. Unpack the P-737 with care.
2. Compare the contents with the items listed in the contract and the packing list.
3. Inspect the contents for signs of damage. If parts are missing or you notice signs of damage, contact PI immediately.
4. Remove the transport lock as described in “Removing the Transport Lock” (p. 2).
5. Keep all packaging materials in case the product needs to be returned.

Removing the Transport Lock

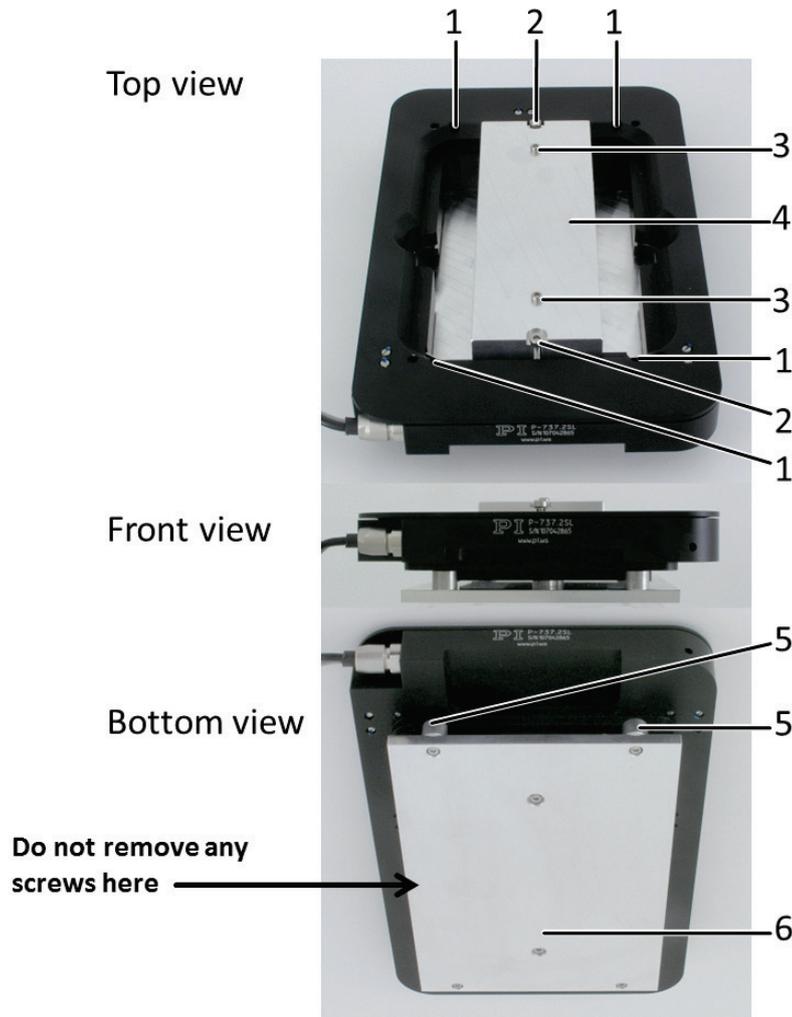


Figure 1: P-737 with transport lock

- 1 M2.5x8 socket head screw (4×)
- 2 M3x16 socket head screw (2×)
- 3 M2.5x20 socket head screw (2×)
- 4 Fixing plate
- 5 Spacer (2 of 6 visible here)
- 6 Base plate

Tools and accessories

- Suitable screwdrivers for M2.5 and M3 socket head screws

Removing the transport lock

1. Remove the fixing plate (4):
To do this, unscrew and remove the two M3x16 screws (2) and the two M2.5x20 screws (3).
2. Remove the base plate (6):
To do this, unscrew and remove the four M2.5x8 screws (1) from the top side.
It is not necessary to remove the six spacers from the base plate.
3. Keep the transport lock in case the product needs to be transported again later.