

MP 80E User Manual  
M-413 Linear Translation Stage  
Release: 1.0.1 Date: 25.04.2007



This document describes the following product(s):

- M-413.xDG  
Linear Translation Stage, DC-Motor
- M-413.xPD  
Linear Translation Stage, ActiveDrive™
- M-413.x2S  
Linear Translation Stage, Stepper Motor



Physik Instrumente (PI) GmbH & Co. KG is the owner of the following company names and trademarks:  
PI®, ActiveDrive™, Mercury™, Mercury™ Step

The following designations are protected company names or registered trademarks of third parties:  
Windows

Copyright 1999–2007 by Physik Instrumente (PI) GmbH & Co. KG, Karlsruhe, Germany.  
The text, photographs and drawings in this manual enjoy copyright protection. With regard thereto, Physik Instrumente (PI) GmbH & Co. KG reserves all rights. Use of said text, photographs and drawings is permitted only in part and only upon citation of the source.

First printing 25.04.2007  
Document Number MP 80E BSc, Release 1.0.1  
M-413\_User\_MP80E\_101.doc

Subject to change without notice. This manual is superseded by any new release. The newest release is available for download at [www.pi.ws](http://www.pi.ws) (<http://www.pi.ws>)

## About This Document

## Users of This Manual

This manual is designed to help the reader to install and operate the M-413 Linear Translation Stage. It assumes that the reader has a fundamental understanding of basic servo systems, as well as motion control concepts and applicable safety procedures.

The manual describes the physical specifications and dimensions of the M-413 Linear Translation Stage as well as the procedures which are required to put the associated motion system into operation.

## Conventions

The notes and symbols used in this manual have the following meanings:



### WARNING

Calls attention to a procedure, practice or condition which, if not correctly performed or adhered to, could result in injury or death.



---

### CAUTION

Calls attention to a procedure, practice, or condition which, if not correctly performed or adhered to, could result in damage to equipment.

---

---

### NOTE

Provides additional information or application hints.

---

The motion controller and the software tools which might be mentioned in this documentation are described in their own manuals. All documents are available as PDF files via download from the PI Website ([www.pi.ws](http://www.pi.ws) (<http://www.pi.ws>)) or on a product CD. For updated releases contact your Physik Instrumente sales engineer or write [info@pi.ws](mailto:info@pi.ws) (<mailto:info@pi.ws>).

# Contents

1	Introduction	3
1.1	Product Description.....	3
1.2	Prescribed Use.....	4
1.3	Safety Instructions.....	5
1.4	Model Survey .....	6
1.5	Unpacking .....	6
1.6	Additional Components .....	7
2	Start-Up	8
2.1	Mechanical Mounting .....	8
2.2	Moving the Stage .....	10
3	Operational Considerations	12
3.1	PWM Amplifiers.....	12
3.2	Limit Switches .....	12
3.3	Reference Switch .....	13
3.4	Rotary Encoder .....	13
3.5	Motor Controllers.....	13
3.6	Host PC.....	13
4	Maintenance	15
5	Troubleshooting	16
5.1	Diagnosis .....	16
5.2	Customer Service.....	18
6	Technical Data	19
6.1	Specifications .....	19
6.2	Range and Transmission Ratio.....	21
6.3	Dimensions .....	21
6.3.1	M-413.xxx .....	22
6.3.2	Accessory Adapter Plates .....	23
6.4	Pin Assignments .....	26
6.4.1	Analog DC-Motor Versions.....	26
6.4.2	DC-Motor ActiveDrive™ (PWM).....	27

6.4.3	2-Phase Stepper Motor .....	27
7	Appendix	28
7.1	Related Products.....	28
7.2	Motor Controller Details .....	29
7.2.1	Stepper motor controllers .....	29
7.2.2	DC Motor Controllers .....	29
8	Glossary of Terms	33
9	Index	37

# 1 Introduction

## 1.1 Product Description



Figure 1: M-413 linear stage versions (from left): the M-413.3PD, M-413.2PD and M-413.1PD provide travel ranges from 100 to 300 mm.

The optimum choice of high-value, low-cost components makes the M-403 family of linear stages very attractive and economical. These stages are leadscrew-driven and provide a minimum incremental motion to 0.2  $\mu\text{m}$ . The precision-machined, high-density, stress-relieved aluminum base provides exceptional stability with minimum weight.

M-413 are available in 3 lengths providing travel ranges from 100 to 300 mm.

### High-Load and Maintenance-Free

All models are equipped with high-precision linear guiding rails and recirculating ball bearings. The recirculating ball bearings are maintenance free and immune to cage migration. The choice of components and careful mounting guarantees high load capacity, longer lifetime and high guiding accuracy.

The stages can carry up to 50 kg and push/pull up to 50 N.

### Three Motor Drive Options

M-413.xPD versions feature the high-performance ActiveDrive™ system. The ActiveDrive™ design, developed by PI, features a high-efficiency PWM (pulse width modulation) servo-amplifier mounted side-by-side with the DC

motor and offers several advantages:

- Increased efficiency, by eliminating power losses between the amplifier and motor
- Reduced cost of ownership and improved reliability, because no external driver is required
- Elimination of PWM amplifier noise radiation, by mounting the amplifier and motor together in a single, electrically shielded case

M-413.xDG models are equipped with a DC motor and a shaft-mounted optical encoder, providing a minimum incremental motion of 0.2  $\mu\text{m}$ .

M-413.x2S models feature a direct-drive, 2-phase stepper motor, providing very smooth operation and a resolution of 0.1  $\mu\text{m}$  per step.

#### Limit and Reference Switches

For the protection of your equipment, non-contact Hall-effect limit and reference switches are installed. The direction-sensing reference switch supports advanced automation applications with high precision.

The M-403 / M-413 and M-404 / M-414 series of linear stages form a modular system. The M-403 is the basic family, providing travel ranges from 25 to 200 mm. M-413 is designed for higher loads with travel ranges from 100 to 300 mm. The M-404 and M-414 stages have the same travel ranges and load capacities, but offer higher precision and more speed..

---

## 1.2 Prescribed Use

Correct operation of the M-413 positioners is only possible in combination with a suitable controller/driver (ordered separately) and software. The controlling device must be able to read out and process the signals from reference and limit switches and the incremental position encoder to ensure proper performance of the servo-control system.

Based on their design and realization, the M-413 stages are intended for single-axis positioning, adjusting and shifting of loads at various velocities. M-413 positioners can be mounted horizontally or vertically. To achieve the specified guiding accuracy, the positioners have to be mounted on a flat surface to avoid torsion of the basic profile.

The positioners may only be used for applications suitably in accordance with the device specifications.

Respect the safety instructions given in this User Manual. The operator is responsible for the correct installation and operation of the M-413.

The verification of the technical specifications by the manufacturer does not imply the validation of complete applications. In fact the operator is responsible for the process validation and the appropriate releases.

The M-413 meets the specifications as defined by DIN EN 61010 for safe operation under normal ambient conditions. See the specifications table for details.

---

## 1.3 Safety Instructions

---

### CAUTION

Read this before operating the equipment covered in this manual. Always keep the User Manual safe and close to the described device. In case of loss or damage to the instructions, please order a new copy from your PI distributor or download one from [www.pi.ws](http://www.pi.ws) (<http://www.pi.ws>). Also keep and add all further information (e.g. extended instructions or Technical Notes) to the User Manual.

---



Read carefully the User Manuals of all other components involved such as controllers or software.

Read the following before operating the equipment covered in this manual:

### WARNING—CRUSH HAZARD

Never put anything where the moving platform or any connected object could possibly trap it! The moving platform can accelerate rapidly and generate high forces which can cause damage or injury.



---

### CAUTION

Be aware that the stage could start an unexpected move when first connected to the motor controller.  
Be aware that failure of the motor controller may crash the stage against a hard stop.



Crashes can cause irreparable damage. Do not disable limit switches in software. Test limit switch operation at low speeds. Stop system if necessary. Do not let it run against a hard stop.

---

## 1.4 Model Survey

A total of 9 M-413 standard versions is available. They differ in terms of drive types and travel ranges. The drive type and the travel range are encoded in the model number.

The drive section options are shown in the table below. The “x” in the model number stands for the travel range code. The major difference between the various models lies in the drive section, as shown in the table below.

Models	M-413.xDG	M-413.xPD	M-413.x2S
Drive type	DC-motor, gearhead, 3.6 W	DC-motor, ActiveDrive™, 42 W	2-phase stepper motor
Position encoder	Yes	Yes	No
Limit switch	Yes	Yes	Yes
Direction-sensing reference switch	Yes	Yes	Yes

The travel range differences and their encoding in the model number are as follows (here the “xx” in the model number stands for the drive-type code):

Models	M-413.1xx	M-413.2xx	M-413.3xx
Travel range in mm	100	200	300

## 1.5 Unpacking

Unpack the M-413 Linear Translation Stage with care. Compare the contents against the items covered by the contract and against the packing list.

The following components are included:

- M-413 Linear Translation Stage
- C-815.38 motor cable for connecting to motor controller
- Screw set with 6 M4 x 30 and 4 M6 x 2 screws for mounting

- User manual for M-413 (this document) in printed form
- M-413.xPD only: M-500.PS power supply with line cord

Inspect the contents for signs of damage. If parts are missing or you notice signs of damage, contact PI immediately.

Save all packing materials in case the product need be shipped again.

---

## 1.6 Additional Components

Depending on your application, you might need the following accessories for use with the M-413 Linear Positioning Stage :

- M-413.AP1: Adapter plate for mounting stages with PI standard hole pattern on M-413 stages,
- M-413.AP2: Z-axis adapter plate for vertical mount of M-413 stages on M-413 stages,
- M-403.AP3: Universal adapter plate for mounting M-413 stages on stages with PI standard hole pattern.

Contact your PI sales engineer or write [info@pi.ws](mailto:info@pi.ws) if you need additional components.

## 2 Start-Up

### 2.1 Mechanical Mounting

#### NOTE

If the counter-sunk mounting holes in the stage base are covered by the platform, you must first move the platform to a position which allows access. Read "Moving the Stage (p. 10)" for instructions.

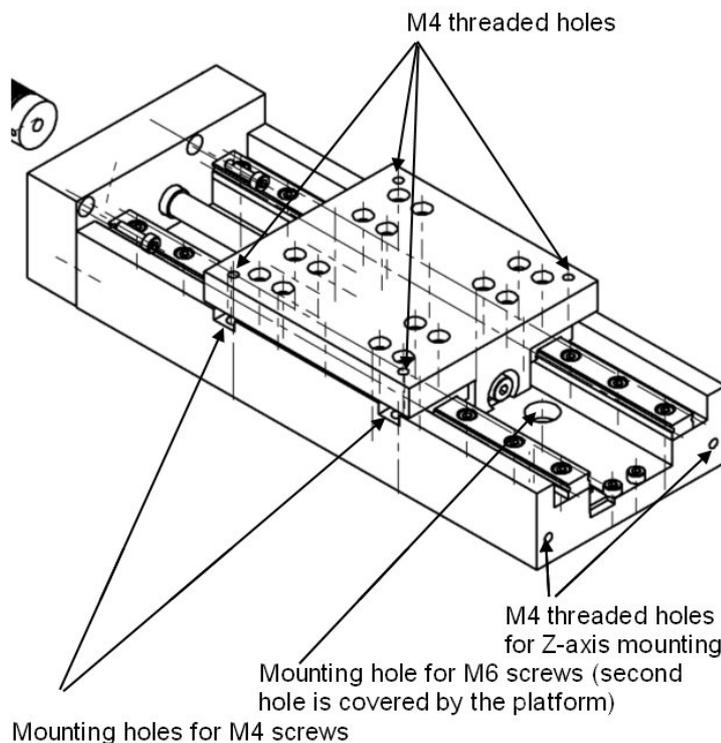



Figure 2: M-ounting hole locations on stage, identical for all models

To mount an M-413 stage on a base plane, use the two holes for M6 screws or the four counter sunk holes for M4 screws in the stage base and the corresponding mounting screws which come with the stage. The arrangement of the M6 holes complies with the hole pattern of standard honeycomb tables.

You have best access to the M6 mounting holes, when the platform is at its reference position (for the referencing procedure see the User Manual of the motor controller used).

You can assemble stacked XY- or XYZ-systems with M-413 stages. If you want to combine different models of the M-403 / M-404 / M-413 / M-414 series, the M-403.AP3 adapter plate might be required. For mounting, use the M4x30 screws which come with the stages:

- For an XY-system, use the four counter-sunk holes for M4 in the base of the upper M-413 stage and the four M4 threaded holes in the moving platform of the lower M-413 stage (see Fig. above).
- For Z-axis mounting, the M-413.AP2 Z-axis adapter plate is required. First mount this plate on the front of the Z-axis stage using the two M4 threaded holes in the stage base (see Fig. above). Then affix the adapter plate to the moving platform of the lower M-413 stage.

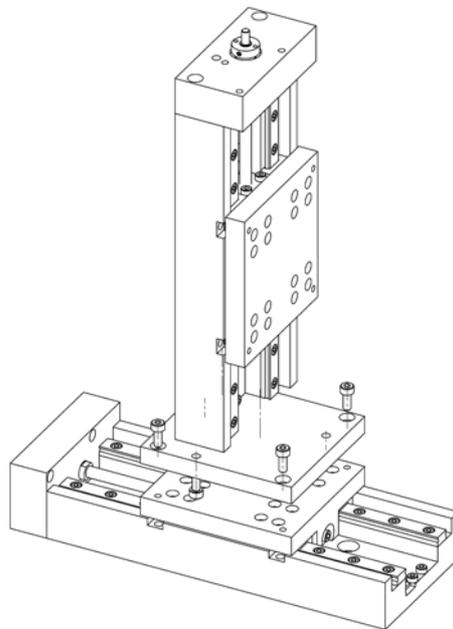

*Figure 3: Z-axis mounting*

To mount a stage with PI Standard Hole Pattern on an M-413 stage, the M-413.AP1 Adapter plate is required.

To mount a M-413 stage upon a stage with PI standard hole pattern, the M-403.AP3 adapter plate is required.

## 2.2 Moving the Stage



### WARNING—CRUSH HAZARD

Never put anything where the moving platform or any connected object could possibly trap it! The moving platform can accelerate rapidly and generate high forces which can cause damage or injury.

### NOTE

The positive direction of motion is away from the end where the motor is mounted.

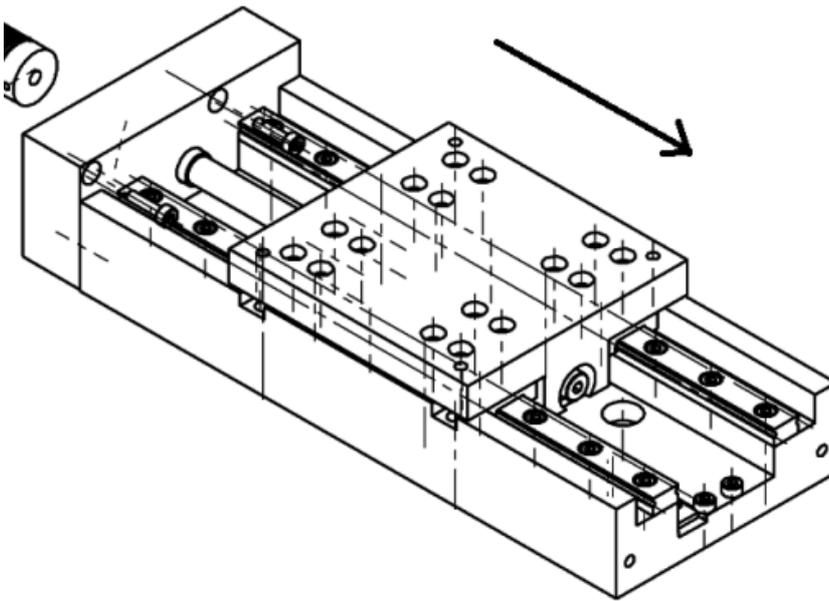


Figure 4: M-4xx stage unit, arrow indicates positive direction of motion

Prior to operation, you should read the documentation for your motor controller. If using any PWM stages read the section about PWM amplifiers (p. 12) carefully.

To operate a stage, proceed as follows:

- 1 Install/connect the controller following the instructions in the controller manual. If you are going to use a host computer, install the

host software in the host computer. This procedure is described in the controller User Manual and/or associated software manuals.

- 2 Connect the stage to the controller using the connecting cable (part number C-815.38), which comes with the stage. With multi-axis controllers, be sure to note the axis designation of the connection selected.
- 3 M-413 versions with ActiveDrive™ (PWM) only: Connect the stage to the M-500.PS power supply, which comes with the stage and connect the power supply to line power (wide range). To activate the M-500.PS power supply switch the "-/o"-button to position "-".
- 4 Command a few test moves to make sure everything is working properly.

For best performance and system compatibility we recommend using PI controllers.

Most PI controllers come with a graphic-user-interface-software which is easy to setup and allows quick installation and testing.

---

## NOTE

Most standard controllers and / or software from PI use a DAT file which contains information on all standard micropositioners and installs automatically with the host software. This file, called pistages.dat, also includes an initial set of PID servo-control parameters for each stage type. If your system does not use pistages.dat, possible PID parameters are included in this manual.

---

## 3 Operational Considerations

### 3.1 PWM Amplifiers

The M-413.xPD stages feature a DC servo-motor with a high-efficiency PWM amplifier mounted side by side. This ActiveDrive™ system provides maximum dynamic performance. An external plug-in power supply (M-500.PS) is provided to supply the built-in amplifier. This architecture allows high torque and high velocities while loading the motor controller with control signals only.

### 3.2 Limit Switches

All M-413 positioners are equipped with non-contact, Hall-effect limit switches (with TTL drivers). Each limit switch sends an overtravel signal on its own dedicated line to the controller. It is the controller that is then responsible for stopping the motion. If it does not do so in time, the positioner will run into the hard stop. On stepper-motor versions, limit switch outputs are active low, on other versions, active high.



#### CAUTION

Crashes can cause irreparable damage. Do not disable limit switches in software. Test limit switch operation at low speeds. Stop system if necessary. Do not let it run against a hard stop.

#### Limit Switch Specifications:

Type:	magnetic (Hall-effect) sensors
Supply.voltage	+5 V / GND, supplied by the motor controller through the motor connector.
Signal output:	TTL level
Sink / Source capab.	20 mA at 18°C
Signal logic (stepper motors)	active low; normal operation: high, limit sensor reached: low
Signal logic (DC-motors):	active-high; normal motor operation: low, limit sensor reached: high

---

### 3.3 Reference Switch

All stage versions are equipped with a direction-sensing Hall-effect reference switch, which is located at about the midpoint of the travel range. This sensor provides a TTL signal indicating whether the positioner is to the positive or negative side of a fixed point. The rising or falling edge of this signal can be used to indicate a known reference position within 0.5  $\mu\text{m}$  accuracy (depending on the controller). The difference in the reference point when approached from the positive or the negative side is about 0.2 mm to 0.4 mm. See the controller User Manual and/or associated software manuals for the commands which make use of the reference signal.

---

### 3.4 Rotary Encoder

M-413.xDG and M-413.xPD are equipped with DC motors with a shaft-mounted position encoder. It provides highly accurate signals whenever the position changes by a known, small amount. By monitoring the encoder pulses, called *counts*, the controller can keep track of the relative motion of the stage. To discover the absolute position, it is necessary to drive the stage to a limit or reference switch. Most controllers have commands to automate this procedure, which is called *referencing*.

---

### 3.5 Motor Controllers

M-413-series micropositioners can be connected to the same motor controllers as other micropositioners from PI. DC motor controllers from PI can control both analog and ActiveDrive™ (PWM) motor versions. The required operating mode is set automatically—depending on the controller model, either directly when the controller is powered up, or when the user chooses the connected positioner type in the control software. If networking several positioners, you have to keep in mind that stepper motor devices can only be networked with other stepper-motor devices. This manual includes a list of suitable controllers.

---

### 3.6 Host PC

Most controllers from PI are in turn controlled by a host PC provided by the user. The controller is either installed directly in the PC (e.g. C-843 PCI board) or connected to it over a communications interface. See the controller User Manual for details.



PI provides software and/or drivers that run on the host PC to control the system. Typically, there will be a control program with a graphic user interface for testing and simple operation, and, in addition, DLL, COM and/or LabView drivers for users who wish to use custom software. If the controller is a PC ISA or PCI card, there will also be hardware drivers to install. See the controller user and software manuals for installation details.

## 4 Maintenance

When operated in a clean environment, no maintenance is required. If the stages are operated in extremely dusty environments, we recommend cleaning and greasing the lead screw/ball screw and linear bearings from time to time. The time interval depends of the degree of contamination and can vary from 100 to 800 operating hours.

# 5 Troubleshooting

---

## 5.1 Diagnosis

Positioner does not move.

Cables not connected properly:

- Check the connecting cables.

For positioners with ActiveDrive™ only:

Check if power supply is connected:

- Verify that the included MS-500.PS power supply is connected and turned on.

Positioner or positioner cable is defective:

- Replace positioner with a working positioner of the same type to test a new combination of controller and positioner (positioners of a different type may malfunction due to unsuitable parameter settings on the controller).

Test cable and / or replace with a working cable of the same type.

Unusually high friction in the spindle-nut-system:

- Clean the spindle and add new lubricant.

Wrong command or wrong syntax:

- Check the error code (with PI-GCS command set, use ERR?; see controller and software manuals for the error code explanations).

Positioner is not referenced:

- In the GCS command set, some closed-loop motion commands are not executed if the mechanics has not been referenced—(ERR? replies "5", see controller and software manuals). Reference the axis as described in the controller and user manuals.

Wrong axis commanded:

- Check if commanded axis is that of the desired positioner.

Incorrect configuration:

- Check for anomalous parameter settings on the controller.

Positioner starts to move but stops after a short time.

Unusually high friction in the spindle-nut-system:

- Clean the spindle and add new lubricant.

Positioner does not reach position with expected accuracy.

Motion parameters are not optimized:

- Adjust the motion parameters according to the values given in this document.

Positioner runs jerkily and not smooth enough.

Motion parameters are not optimized:

- Adjust the motion parameters according to the values listed in this document.

Positioner did not stop in time and ran into a hard stop.

The controller has not stopped the positioner in time:

- Too high velocity. See chapter "Limit Switches".
- Turn off the motor. If possible, you may push the positioner out of the end zone manually.
- The end of the leadscrew has an Allan wrench socket so it can be cranked manually.

---

## 5.2 Customer Service

Still having problems? Call your PI representative or write to [info@pi.ws](mailto:info@pi.ws); please have the following information about your system ready:

- Product codes and serial numbers of all products in the system
- Current firmware version of the controller (if present)
- Software version of drivers and / or host software (if present)
- Operating system on host PC (if present)

# 6 Technical Data

## 6.1 Specifications

Models	PD (ActiveDrive™) versions M-413.1PD/M-413.2PD M-413.3PD	DG (gearhead) versions M-413.1DG/M-413.2DG M-413.3DG	2S (stepper) versions M-413.12S/M-413.22S M-413.32S	Units	Tolerance
Active Axes	X	X	X		
Motion and Positioning					
Travel range	100 / 200 300	100 / 200 300	100 / 200 300	mm	
Integrated sensor	Rotary encoder	Rotary encoder	—		
Sensor resolution	4000	2000	—	cts./rev.	
Design resolution	0.25	0.018	0.16	µm	typ.
Min. incremental motion	0.25	0.2	0.2	µm	typ.
Backlash	6	10	6	µm	typ.
Unidirectional repeatability	1	1	1	µm	typ.
Pitch	300	300	300	µrad	typ. over 100 mm
Yaw	300	300	300	µrad	typ. over 100 mm
Max. velocity	10 <sup>#</sup>	2.5	3	mm/s	
Origin repeatability	1	1	1	µm	typ.
Mechanical properties					
Spindle pitch	1	1	1	mm/U	
Gear ratio	—	28.444	—		

Motor resolution	—	—	6400*	steps/rev.	
Stiffness in motion direction	6000	6000	6000	N/mm	± 20%
Load Capacity**	500	500	500	N	max.
Push / pull force	50	50	50	N	max.
Lateral Force	200	200	200	N	max.
Drive Properties					
Drive type	DC Motor, ActiveDrive™	DC motor, gearhead	2-phase stepper motor*		
Operating voltage	24 (PWM)	0-12 differential, polarity determines direction	24	V	
Electrical power	70	3.6	4.8	W	nominal
Miscellaneous					
Operating temperature range	-20 bis +65	-20 bis +65	-20 bis +65	°C	
Material	Al (black anodized)	Al (black anodized)	Al (black anodized)		
Mass	4.4 / 5.4 6.6	4.2 / 5.2 6.4	4.4 / 5.4 6.6	kg	± 5%
Cable length	3	3	3	m	± 10 mm
Connector	Sub-D 15 (m)	Sub-D 15 (m)	Sub-D 15 (m)		
Recommended controller/driver	C-862 single-axis C-843 PCI board, for up to 4 axes	C-862 single-axis C-843 PCI board, for up to 4 axes	C-663 single-axis		

\* \*\* 2-phase stepper motor, 24 V chopper voltage, max. 0.8 A/phase, 400 full steps/rev., motor resolution with C-663 stepper motor controller

\*\* Load: Max. normal load capacity: Centered, vertical load (horizontal installation of the stage).

# Max. recommended velocity




## 6.2 Range and Transmission Ratio

	Range [mm]	Range [cts]	Transmission Ratio
M-413.1PD	100	400,000	4 cts/μm
M-413.2PD	200	800,000	
M-413.3PD	300	1,200,000	
M-413.1DG	100	5,688,888	56.88889 cts/μm
M-413.2DG	200	11,377,776	
M-413.3DG	300	17,066,664	

## 6.3 Dimensions

Dimensions in mm, decimal places separated by commas in drawings

**6.3.1 M-413.xxx**

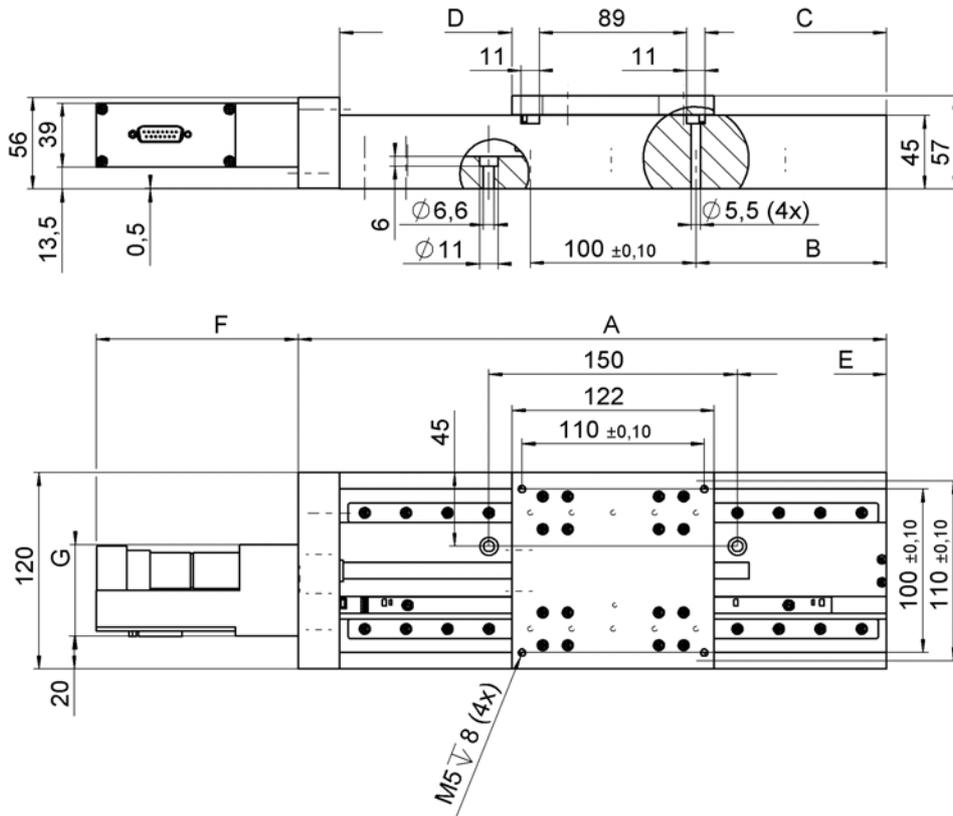

	M-413.1xx	M-413.2xx	M-413.3xx
A	255	355	455
B	65	115	165
C	59,5	109,5	159,5
D	54 ( $\pm 50$ )	104 ( $\pm 100$ )	154 ( $\pm 150$ )
E	40	90	140
	M-413.xPD	M-413.xDG	M-413.x2S
F	120	122	90
G	59,5	56	59,5

### 6.3.2 Accessory Adapter Plates

M-413.AP1

Adapter plate for mounting stages with PI Standard Hole Pattern on M-413 stages

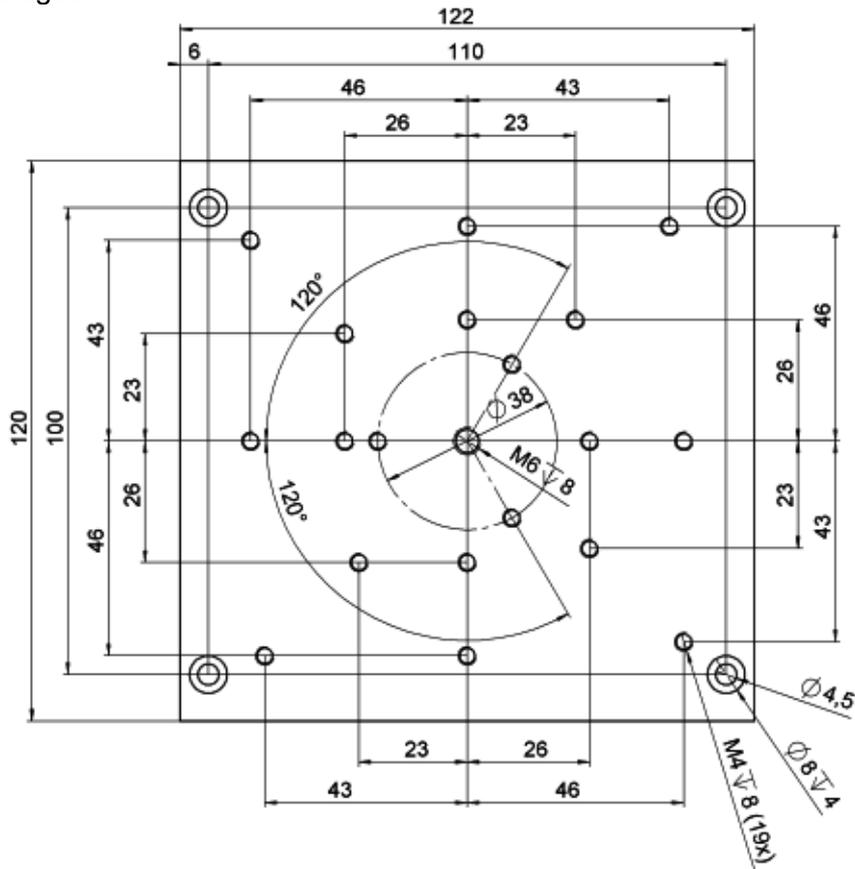


Figure 5: M-413.AP1 adapter plate with PI standard hole pattern

M-413.AP2

Z-axis adapter plate for vertical mount of M-413 stages on M-413 stages

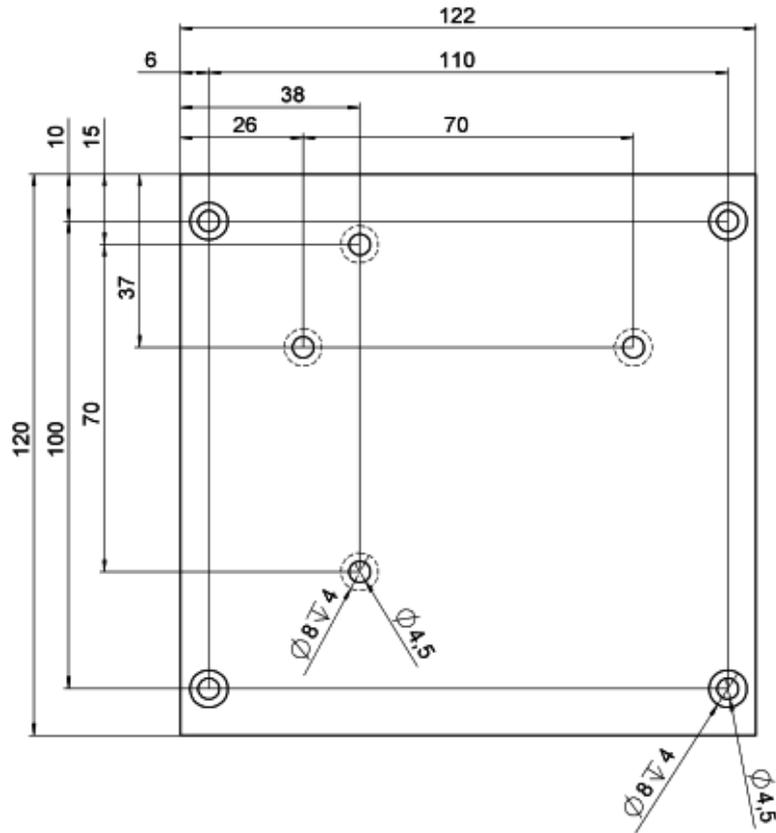


Figure 6: M-413.AP2 Z-axis adapter plate

M-403.AP3

Universal adapter plate for mounting M-413 stages on stages with PI standard hole pattern

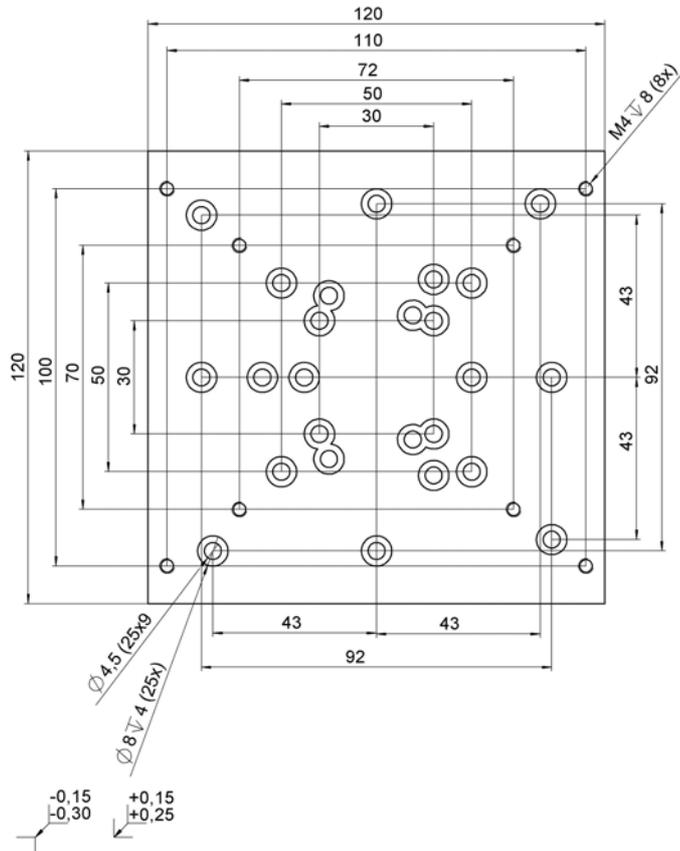


Figure 7: Universal adapter plate



## 6.4 Pin Assignments

### 6.4.1 Analog DC-Motor Versions

Connector type on positioner: D-Sub15m

Pin#	Signal	Direction	Function
1	n.c.		not connected
9	MOT (-)	input	motor connection (-)*
2	MOT (+)	input	motor connection (+)*
10	PGND	input	ground (power)
3	n.c.		not connected
11	n.c.		not connected
4	+5V	input	+5 V input for encoder and logic
12	NLIM	output	negative limit signal (active high), TTL
5	PLIM	output	positive limit signal (active high), TTL
13	REFS	output	position reference signal, TTL
6	GND		ground (logic)
14	A (+)	output	encoder signal A, TTL
7	A (-)	output	encoder signal A-dash, TTL
15	B (+)	output	encoder signal B, TTL
8	B (-)	output	encoder signal B-dash, TTL

\* Polarity determines direction of motion: if MOT+ > MOT- motion is in the positive direction, i.e. away from the end where the motor is mounted.



### 6.4.2 DC-Motor ActiveDrive™ (PWM)

Connector type on positioner: D-Sub15m

Pin#	Signal	Direction	Function
1	ENABLE	input	Enable signal for power amplifier and motor brake, (+5 V to +12 V)
9	n.c.		Not connected
2	n.c.		Not connected
10	PGND	input	Ground (Power)
3	MAGN	input	PWM signal magnitude (speed)
11	SIGN	input	PWM signal sign (direction)
4	+5V	input	+5 V input for encoder and logic
12	NLIM	output	Negative limit signal (active high), TTL
5	PLIM	output	Positive limit signal (active high), TTL
13	REFS	output	Position reference signal, TTL
6	GND		Ground (logic)
14	A(+)	output	Encoder signal A, TTL
7	A(-)	output	Encoder signal A-dash, TTL
15	B(+)	output	Encoder signal B, TTL
8	B(-)	output	Encoder signal B-dash, TTL

### 6.4.3 2-Phase Stepper Motor

Connector type on positioner: D-Sub15m

Pin #	Signal	Direction	Function
1	phase 1a	input	motor winding 1a
9	phase 1b	input	motor winding 1b
2	phase 2a	input	motor winding 2a
10	phase 2b	input	motor winding 2b
3	n.c.		not connected
11	n.c.		not connected
4	n.c.		not connected
12	n.c.		not connected
5	n.c.		not connected
13	n.c.		not connected
6	+5V	input	+5V supply from controller
14	PLIM	output	positive limit signal (active low), TTL
7	GND		ground (power)
15	REFS	output	reference signal, TTL
8	NLIM	output	negative limit signal (active low), TTL

## 7 Appendix

### 7.1 Related Products

With the M-403/M-413 linear stage series PI offers a flexible and cost-effective product line for cost-sensitive applications. The flexibility in travel range (25 to 300 mm) and motorization (stepper motors, gearbox and ActiveDrive™ direct-drive DC motors) provides a broad range of solutions for standard micropositioning tasks.

For more-demanding applications, PI offers the M-404/M-414 stage series. Their outer dimensions and travel ranges are identical to the M-403/M-413, but due to the high-precision components used, they provide improved features such as:

- More accurate guiding system by additional polishing of the bearing surface,
- DC-motor with low-backlash gearhead,
- Higher velocities and duty cycles due to the low friction of the ball screw,
- Higher mechanical stability due to 2 polished mounting steps for low surface contact.

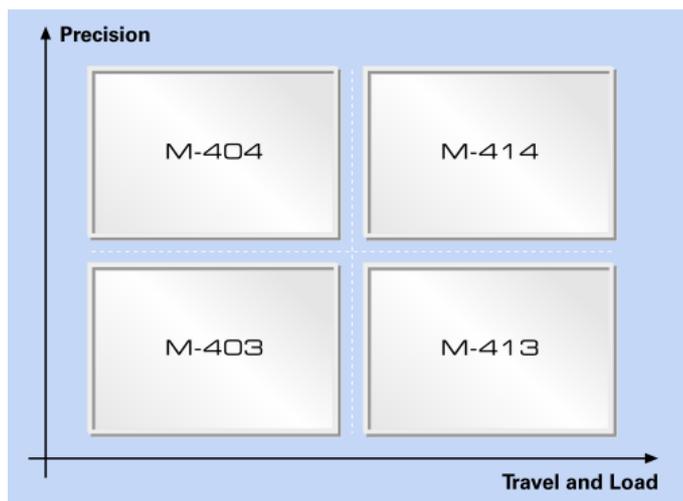


Figure 8: M-403 to M-414 product portfolio. The M-403/M-413 series provide especially cost-effective micropositioning solutions.

The smaller M-403 and M-404 provide a profile width of 80 mm and travel ranges of 25, 50, 100, 150 and 200 mm.  
 The heavy-duty M-413 and M-414 provide a profile width of 120 mm and travel ranges of 100, 200 and 300 mm.

## 7.2 Motor Controller Details

Each motorized stage in the motion system must be connected to a motion controller. The controller is either preprogrammed for stand-alone operation or networked with or installed in a PC: i.e. controller setup and/or operation are effectuated with software.

### 7.2.1 Stepper motor controllers

The M-413 .x2S (stepper motor version) can be controlled with the C-663 Mercury™ Step stepper motor controller. PI stepper motors operate in open-loop mode.

Drive type	Stepper Motor
Controller	C-663 Mercury™ Step
Axes per controller	1
Host PC interface	RS-232 bus or daisy chain
Multiple controllers on same host PC	yes, same port

### 7.2.2 DC Motor Controllers

M-413 DC-motor versions can be used with C-843 motor controller card or C-862 Mercury™ controller.

Drive type	DC Motor	
Controller	C-862 Mercury™	C-843 motor controller card
Axes per controller	1	2 or 4
Host PC interface	RS-232 bus or daisy chain	Internal (PCI bus)
Multiple controllers on same host PC	yes, same or separate ports	yes, separate slots

#### NOTE

Most standard controllers and / or software from PI use a DAT file which contains information on all standard micropositioners and installs

automatically with the host software. This file, called `pistages.dat`, also includes an initial set of PID servo-control parameters for each stage type. If your system does not use `pistages.dat`, possible PID parameters are included in this manual.

---

### Parameter settings for C-862 Mercury™ DC-Motor Controller

The C-862 Mercury™ Controller is a compact palm-size, single-axis controller. C-862s are daisy-chain networkable and a multi-axis system can be controlled from a single PC RS-232 interface.

The Windows operating program for the C-862 Mercury™ controller allows choice of micropositioners as a start-up option for operation. The C-862 Mercury™ has the correct operating mode (analog or PWM) automatically enabled with no extra command required.

### Appropriate PID-Parameters for M-413.xDG

Parameter	Recommended Value	Minimum Value	Maximum Value
p-term	280	40	450
i-term	35	0	150
d-term	250	0	400
i-limit	2000	0	2000
velocity [cts/s]	112,000	0	143,000
velocity [mm/s]	2.0	0	2.5
acceleration [cts/s <sup>2</sup> ]	1,000,000	5,600	2,000,000
acceleration [mm/s <sup>2</sup> ]	18	0.1	36

### Appropriate PID-Parameters for M-413.xPD

Parameter	Recommended Value	Minimum Value	Maximum Value
p-term	220	40	400
i-term	40	0	150
d-term	250	0	400
i-limit	2000	0	2000
velocity* [cts/s]	35,000	0	40,000
velocity* [mm/s]	8.75	0	10
acceleration [cts/s <sup>2</sup> ]	400,000	5,000	1,200,000
acceleration [mm/s <sup>2</sup> ]	100	1.25	300

\* Higher velocities up to 100,000 cts/s (25 mm/s) are possible, but not recommended for continuous operation due to frictional heat generation in the spindle system. Setting velocity limit to 0 has the obvious effect of preventing all motion with no error indications.



## 8 Glossary of Terms

### *ActiveDrive™ DC Motor*

Some of the advantages of DC-motor drives are good dynamic performance, fast response, high torque at low rpm, low heat dissipation and low vibration. The cost of a high-performance amplifier, however, is generally higher than that for a stepper motor.

The ActiveDrive™ system reduces this cost considerably, by integrating a PWM (pulse width modulation) driver and amplifier in the motor case. This design provides several advantages:

- Increased efficiency, by eliminating power losses between the amplifier and motor
- Reduced cost, more-compact system, and improved reliability, because no external driver and cabling are required
- Elimination of PWM amplifier noise radiation by mounting the amplifier and motor together in a single electrically shielded case

Positioning accuracy is assured with either rotary or optical linear encoders.

### *Backlash*

Position error that appears upon reversing direction. Backlash is caused by play in the drive train components coming after the encoder, such as gearheads or bearings, and by friction in the guiding system. Unlike hysteresis, it can lead to instability in closed-loop setups because it causes a deadband in the servo-loop. Some manufacturers promote controllers with automatic backlash compensation that add the estimated amount of lost motion upon each reversal. This solution is very limited in practice, as backlash is not constant but varies with temperature, deceleration, acceleration, load, leadscrew position, direction, wear, etc.

### *Bidirectional Repeatability*

The accuracy of returning to a position from any position, regardless of direction. Effects such as hysteresis and backlash affect bidirectional repeatability. See also "Unidirectional Repeatability".

### *DC Servo-Motor*

A direct-current motor that is operated in a closed-loop system (servo-loop). Characteristics of DC servo-motors are lack of vibration, smooth running, wide speed range and very good low-speed torque. For optimum performance, a good motor controller with PID (proportional, integral, derivative) algorithm and filter settings is mandatory.

### *Design Resolution*

The theoretical minimum movement that can be made. This is a calculated value based on the drive components (drive screw pitch, gear ratio, motor angular resolution etc.) and does not account for nonlinearities like friction, backlash, etc. Design resolution must not be confused with minimum incremental motion. In systems with high gear ratios or microstepping motors, the design resolution can be in the subnanometer range. In practice, incremental motion of less than 0.1  $\mu\text{m}$  is prevented by guiding system friction (except with air bearings and flexures).

### *Maximum Push/Pull Force*

Active and passive force limit in operating direction, at center of stage. Some stages may be able to generate higher forces at the cost of reduced lifetime.

### *Minimum Incremental Motion*

The minimum motion that can be repeatedly executed for a given input, which is sometimes referred to as practical or operational resolution. Design resolution and practical resolution must be distinguished. Design resolutions of 1 nm or better can be achieved with many motor, gearbox and leadscrew combinations. In practical applications, however, stiction/friction, windup, and elastic deformation limit resolution to fractions of a micron. Minimum incremental motion must be determined by actual measurements.

Repeatable nanometer or sub-nanometer resolution can be provided by solid-state actuators (PZTs) and PZT flexure stages (see the "PZT Flexure NanoPositioners" and "PZT Actuators" sections of the PI Catalog for details) such as those integrated in PI's M-511.HD and M-714 Hybrid Drive stages.

### *Pulse Width Modulation (PWM)*

The PWM mode is a highly effective method of transmitting electrical energy at a variable rate by varying the width of pulses in a train rather than the amplitude of an analog signal.

## *Stepper Motor*

An electric motor providing motion in discrete angular steps, without the need of position sensor or servo-loop. Compared to closed-loop DC motors of the same size, stepper motors provide poorer dynamic performance and dissipate more heat, especially in steady state operation. Their advantages lie in long lifetime and simple control electronics. Most common are 2- and 4-phase designs, which are very inexpensive; 3- and 5-phase motors offer better performance in terms of dynamics, torque, stiffness, resolution and noise generation, but their manufacture is more complicated.

## *Unidirectional Repeatability*

The accuracy of returning to a given position from the same direction. Because unidirectional repeatability is almost unaffected by backlash and hysteresis, it is often considerably better than bidirectional repeatability.





# 9 Index

## 2

2-Phase Stepper Motor • 6

## A

Accessory Adapter Plates • 5  
ActiveDrive™ DC Motor • 7  
Additional Components • 2  
Analog DC-Motor Versions • 5  
Appendix • 6

## B

Backlash • 7  
Bidirectional Repeatability • 7

## C

Customer Service • 2

## D

DC Motor Controllers • 7  
DC Servo-Motor • 7  
DC-Motor ActiveDrive™ (PWM) • 6  
Design Resolution • 7  
Diagnosis • 2  
Dimensions • 4

## H

Host PC • 2

## I

Introduction • i

## L

Limit Switches • 2

## M

M-413.xxx • 4  
Maintenance • 2  
Maximum Push/Pull Force • 7  
Mechanical Mounting • 2  
Minimum Incremental Motion • 7

Model Survey • i  
Motor Controller Details • 6  
Motor Controllers • 2  
Moving the Stage • 2

## O

Operational Considerations • 2

## P

Pin Assignments • 5  
Prescribed Use • i  
Produktbeschreibung • i  
Pulse Width Modulation (PWM) • 7  
PWM Amplifiers • 2

## R

Range and Transmission Ratio • 4  
Reference Switch • 2  
Related Products • 6  
Rotary Encoder • 2

## S

Safety Instructions • i  
Specifications • 2  
Start-Up • 2  
Stepper Motor • 7  
Stepper motor controllers • 6

## T

Technical Data • 2  
Troubleshooting • 2

## U

Unidirectional Repeatability • 7  
Unpacking • 2